F · S · T

EN

DE

CN

ES

FR

CARE INSTRUCTIONS FOR YOUR NEW SPRING SCISSORS

FINE SCIENCE TOOLS

THANK YOU FOR BUYING A PAIR OF HIGH-QUALITY FST® SPRING SCISSORS

Every single one of our Spring Scissors has been tested by our surgical technicians and adapted to your needs at the micrometer level.

The result: extremely fine cutting edges, joints and screws – all of which require the same level of care when it comes to handling and maintenance as they did when they were being produced.

- You should clean, rinse and dry the instruments carefully after use.
- For cleaning, we recommend a soft toothbrush, always brushing towards the tip to remove any visible debris, especially from hinges, joints, and other hard-to-reach areas.
- After every other sterilisation, you should wet the screws, joints and moving parts of the scissors with a drop of our instrument oil FST Item No. 29055-00. Please do not use WD-40 or other common industrial oils for this.

This way, you can avoid corrosion and damage to the fine springs, rongeurs, scissors, needle holders and clamps – and ensure that you can always 100% rely on your instruments.

If you have any other technical questions, please don't hesitate to contact us.

STERILISATION TIPS

STEAM STERILISATION / AUTOCLAVE PROCESS

The autoclave process uses pressurised moist heat for sterilisation. It's important that the steam used comes from sterilised water and is not contaminated.

STERILISATION TIPS

- You can sterilise your instruments individually or as a set.
 - Individual instruments:
Disposable cases or sterilisation bags are ideal for this (FST Item No. 28000-06, 28000-11, 28000-14). Make sure that the bag is sufficiently large – the instrument should be in an open (unlocked) position.
 - Instrument sets:
The instruments should all be open/unlocked here too. If it is necessary to have two layers, the heavy instruments should be at the bottom. Otherwise, the steam won't be able to reach and sterilise the adjoining metallic surfaces. In addition, thermal expansion during autoclaving may result in cracks in the joint areas.
- IMPORTANT:** Please DO NOT mix instruments of different or unknown alloys in the autoclave.

STERILISATION TIPS

- Do not overload the autoclave chamber. Doing so would impede steam penetration, meaning the instruments may not be 100% sterilised.
- Place a towel beneath the pan to catch excess moisture during autoclaving. Make sure that the towel has no washing detergent residue on it, and is pH-neutral when submerged in water.
- At the end of the autoclaving cycle (before the drying cycle), unlock the autoclave door and open it slightly. Then allow the drying cycle to run for the period of time recommended by the autoclave manufacturer.

PLEASE NOTE: If you open the autoclave door fully before the drying cycle, cold air will flow into the chamber – and this may lead to condensation forming on the instruments and cause water spots on the instruments.

STERILISATION TIPS

ETHYLENE OXIDE STERILISATION (ETO STERILISATION)

Some instruments, such as catheters, cannulas, etc., cannot withstand high temperatures or high humidity levels. Ethylene oxide sterilisation (ETO sterilisation) is recommended for these instruments. This is a process composed of four primary variables: gas concentration, humidity, temperature and time.

You can find the appropriate sterilization accessories under the following product numbers in our webshop or catalog:
FST Item No. 28000-06, 28000-11, 28000-14, 20840-01, 20840-02, 20840-03, 20842-03, 20810-01, 20810-02, 20850-00

OXIDATION (RUST)

FST provides high quality, stainless steel surgical instruments that are resistant to oxidation or rust. Each individual instrument is manufactured in strict accordance with European standards for stainless steel medical goods.

You should maintain these instruments with great care in order to extend their service life. Stainless steel is, of course, corrosion-resistant. Nevertheless, improper handling may result in rusting or discolouration – if stainless steel is subjected to physical, thermal or chemical damage, it may lose some of its rust-proof properties.

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