F · S · T

EN

DE

CN

ES

FR

INSTRUCTION MANUAL AUTOCLIP® 9 mm KIT

FINE SCIENCE TOOLS

AUTOCLIP® 9 mm KIT

FST® ITEM NO. 12020-00

LOADING THE APPLIER

1. Compress the spring assembly using your thumb, then flip the spring retainer of the applicator and remove the spring assembly.

2. Insert clips arrow end first.


3. Reinsert the spring assembly and close the spring retainer.

4. Pull out the blue clip rack. If removal is difficult, carefully squeeze the prongs of the rack together while removing it.

THE APPLIER IN USE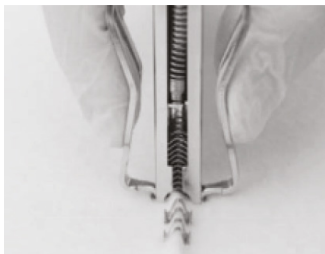

9 mm AutoClip® Wound Clip

1. Hold the applicator as close to its jaws as convenient. The applicator has indentations for correct finger and thumb positioning and should be held at the depressions.
2. Place the applicator jaws on the outer side of the wound to be closed. Make sure the tissue layers are aligned for secure and proper placement.
3. Apply the clip by squeezing the applicator quickly and firmly.
4. After a clip has been applied, allow the jaws to open fully by releasing the pressure.

PRECAUTIONS

Treat any component of the system as potentially bio-hazardous after contact with blood or tissue. Use universal precautions and safe laboratory practices.

STERILIZATION

Use your established procedures to sterilize the AutoClip Wound Clip Applicator and AutoClip Wound Clips.

If a sterilization procedure has not been established, we recommend using wet steam sterilization at 120 – 125 °C for at least 30 minutes in an autoclave to sterilize the device.

REMOVING THE CLIPS

1. Hold the remover at a 45° angle and align the remover tips with the wound lip grooves.
2. Compress the remover to open the wound clips.

NOTE: Unused clips may be recovered from the applicator by reinserting the carrier rack into the remaining clips. Then remove the spring assembly and carrier rack. Clips from more than one rack may be combined to make fully loaded racks of twenty.

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