F · S · T

EN

DE

CN

ES

FR

PRODUCT MANUAL HOT BEAD STERILIZER

FINE SCIENCE TOOLS

HOT BEAD STERILIZER

FST® Item No. 18000-45 – Well Volume: 100 ml

FST® Item No. 18000-50 – Well Volume: 175 ml

INSTRUCTIONS

Fill the sterilizer well with the supplied glass beads to within 3 mm below the rim.

Insert electrical plug into wall socket. DO NOT cover well with lid when unit is hot. Switch unit on. It will take 15 to 20 minutes to reach operating temperature of approximately 240 to 270 °C.

Carefully insert the working part of clean surgical instrument into well after unit has reached operating temperature and leave for 10 to 20 seconds. Do not leave longer as instruments become too hot to handle. Sterilization is on contact with glass beads, so tubes (cannulae) cannot be sterilized effectively.

Unit can be left in the “ON” position all day and is safe for use in laminar flow cabinets.

Only when unit is switched off and cool, cover with lid to reduce dust accumulation. Should unit overheat and trip safety switch, let it cool off completely, then turn it over and push in red button visible inside the unit as viewed through one of the cooling holes.

INSTRUCTIONS

The glass beads should be cleaned routinely and replaced in a dry state. Use a spoon for removal of the glass beads, since otherwise there is a slight chance of beads getting stuck in the rim of the well.

IMPORTANT: This unit should only be used for research purposes. DO NOT use in situations which involve humans or valuable animals. DO NOT USE in surroundings where inflammable gases / substances are involved.

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