F · S · T

EN

DE

CN

ES

FR

INSTRUCTION MANUAL MAGNETIC FIXATOR RETRACTION SYSTEM

FINE SCIENCE TOOLS

MAGNETIC FIXATOR RETRACTION SYSTEM

FST® Item No. 18200-50 – Kit for animals from 200 g up to 5 kg

FST® Item No. 18200-20 – Kit for animals up to 200 g

Included with kit:

Magnetic stand (18200-01, 18200-02) **A**


Base Plate (18200-03, 18200-04) **B**

Wire Handles (18200-05, 18200-06) **C**

Elastomer (18200-07) – not autoclavable; for single use only! **D**

Sharp Retractors (18200-08) **E**


Blunt Retractors (18200-09, 18200-10, 18200-11, 18200-12) **F**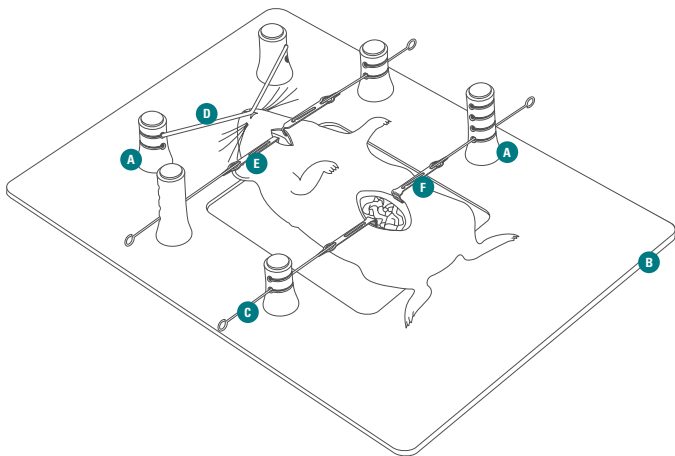

BASE PLATES

Base plates are available in small (FST Item No. 18200-03) and large (FST Item No. 18200-04) sizes. The small base plate is ideal for mouse surgeries, and the large base plate is best suited for rat, guinea pig and small rabbit procedures. Both have an internal window which allows the animal to rest upon either a temperature maintenance system or insulating material.

The base plates are made of ferromagnetic stainless steel and can be treated like other stainless steel trays.

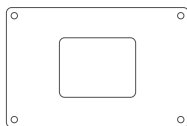
Small (20 x 30 cm)

FST Item No. 18200-03

Large (25 x 35 cm)

FST Item No. 18200-04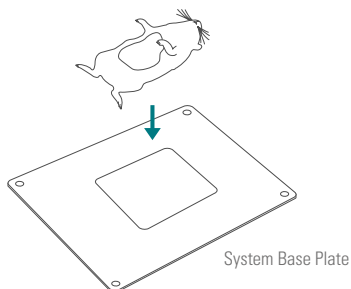


MAGNETIC FIXATORS

Fixators are used to hold the retractors. They contain rare earth magnets that attach anywhere on the base plate. A push button top operates multi-level locking jaws that grip wire or elastomer firmly. They can be adjusted linearly or rotationally with finger-tip release.

The magnetic field is well controlled within the fixator to minimize interference with sensitive instrumentation.

Magnetic fixators come in two sizes. The short fixator (FST Item No. 18200-01) offers two locking slots and is ideal for small animal procedures or where a low profile fixator is desired. The tall fixator (FST Item No. 18200-02) offers four locking slots and is best suited for large animal procedures.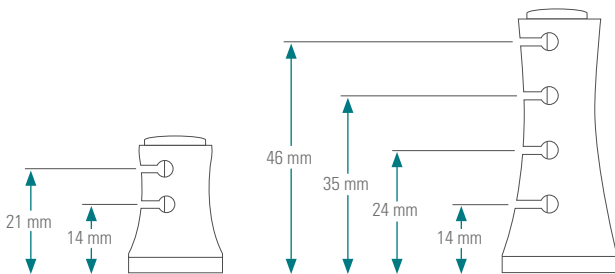


Short

FST Item No. 18200-01


Tall

FST Item No. 18200-02

MAGNETIC FIXATORS
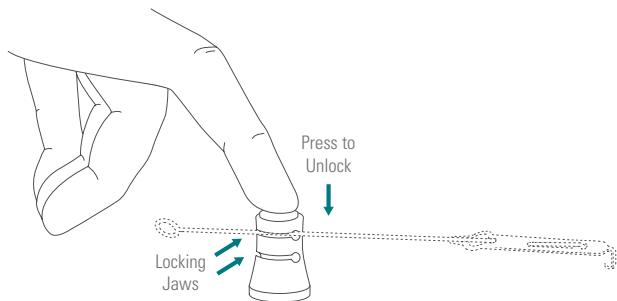


ANATOMY OF THE FIXATOR

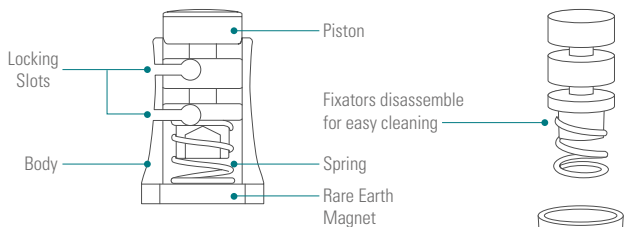CAUTION!




Rare earth magnet in base should not be exposed to temperatures in excess of 148 °C.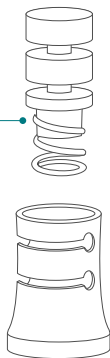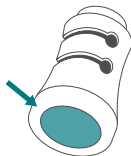

RETRACTOR COMPONENTS


RETRACTOR TIPS

Tip Profile

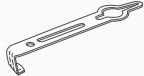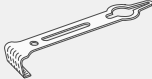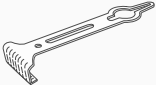
		SHARP use as a skin hook or as a traumatic tissue retractor FST Item No. 18200-08
		BLUNT 1 mm FST Item No. 18200-09
		BLUNT 2.5 mm FST Item No. 18200-10
		BLUNT 5 mm FST Item No. 18200-11
		BLUNT 7.5 mm FST Item No. 18200-12

RETRACTOR COMPONENTS

WIRE HANDLES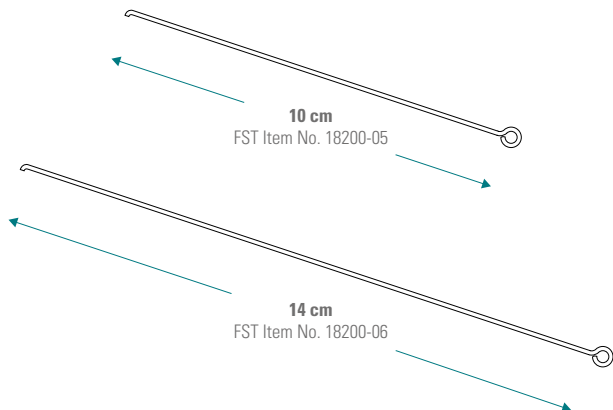

ELASTOMER

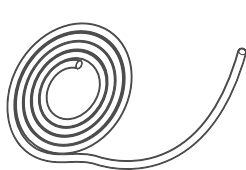FST Item No. 18200-07


CAUTION!

Not autoclavable!


STATIC RETRACTION

Static retraction uses wires to provide precise positioning of the wound margins.

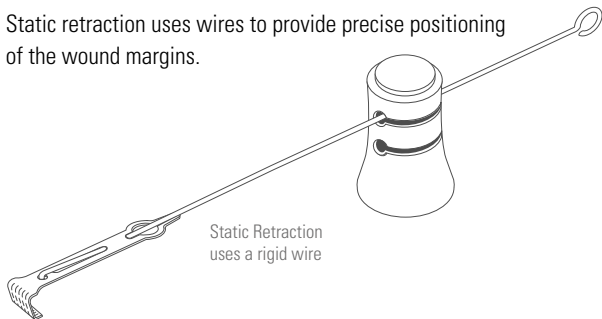
A 10 cm or 14 cm wire handle interlocks with any retractor tip to form a static retractor.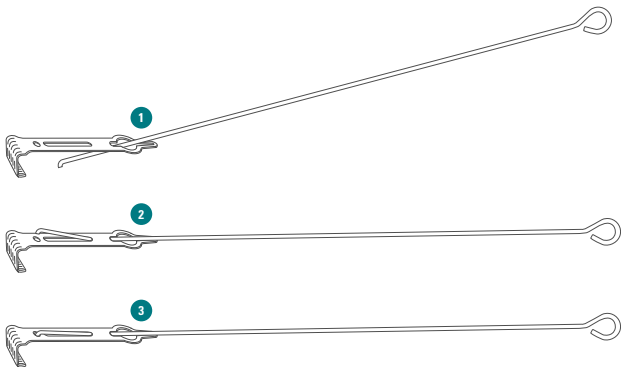


DYNAMIC RETRACTION

Dynamic retraction uses elastomer to provide a constant force on the wound margins.

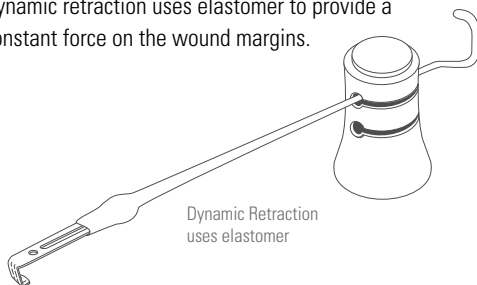
A length of elastomer can be over-sleeved (slid over) any retractor cleat to create permanent dynamic retractors.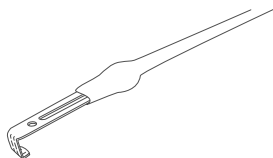


OR

A length of elastomer can be locked into the internal cleat of any retractor.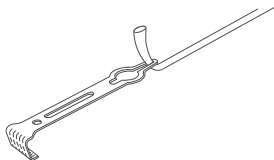


THE SYSTEM IN USE

Either a wire or elastomer can be secured in the jaws of the fixator, and easily adjusted or retensioned at any point during the procedure.

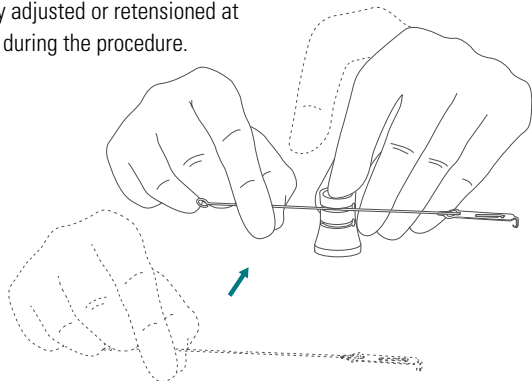
Fixators can also be repositioned at any time during the procedure.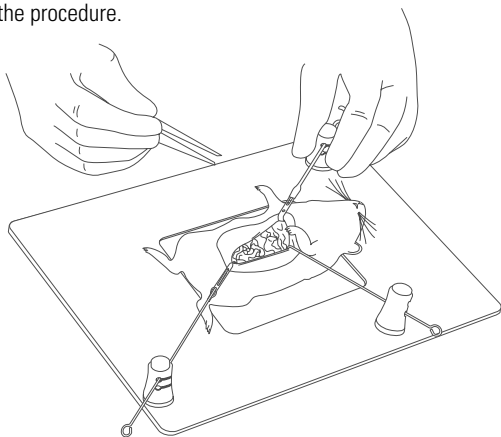

THE SYSTEM IN USE

Other components such as anesthesia inhalation cones can be wrapped in monofilament and positioned using the locking jaws of the fixators. Wires for devices such as temperature probes can be held directly by the fixator jaws.

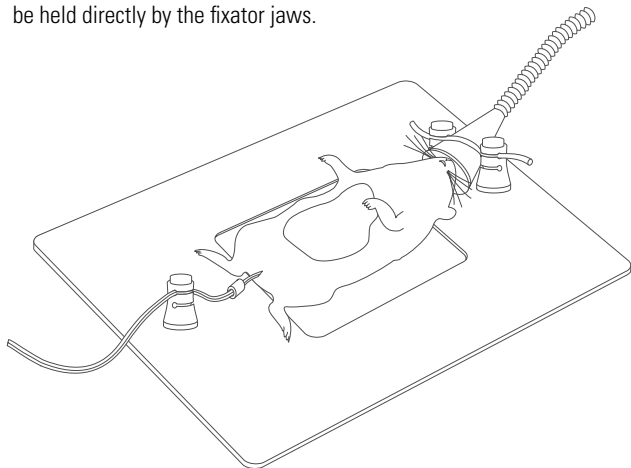AVAILABLE COMPONENTS

FST Item No.	Name
18200-03	Small base plate (20 cm x 30 cm)
18200-04	Large base plate (25 cm x 35 cm)
18200-01	Short fixator (height: 2.5 cm)
18200-02	Tall fixator (height: 5 cm)
18200-05	Wire handles (length: 10 cm; Pkg. of 10)
18200-06	Wire handles (length: 14 cm; Pkg. of 10)
18200-07	Elastomer (2 m roll)
18200-08	Retractor, sharp point (Pkg. of 10)
18200-09	Retractor, blunt (width: 1 mm; Pkg. of 10)
18200-10	Retractor, blunt (width: 2.5 mm; Pkg. of 10)
18200-11	Retractor, blunt (width: 5 mm; Pkg. of 10)
18200-12	Retractor, blunt (width: 7.5 mm; Pkg. of 10)

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