



Hot Bead Sterilizer - 8cm Well



This sterilizer provides a safe, effective and convenient method for sterilizing your surgical instruments prior to surgery or tissue culture. Without using potentially dangerous gases, flames or chemicals, you can ensure that your instruments will be free of pathogens and microbial contaminants. The beads take 15-20 minutes to reach 240°C to 270°C (464°F to 518°F), and then the sterilizer is ready for use. The temperature is thermostatically maintained, and the unit can be left on all day. Instruments require approximately 20 seconds of contact with the beads for complete sterilization. The sterilizer is safe to use even in laminar flow hoods. Instruments not included.

General Info

Product Number	18000-45
Category	Instrument Care > Sterilization
EAN	7612345006789



Find 18000-45 on
finescience.com

Technical Info

Shipping Weight	1770 g	Net Weight	1530 g
Package Size	1 Pc.	Unit of Measure	EA
Minimum Order Quantity	1		
Alloymaterial	Various		
Dimensionstext	14.5 x 12.5 x 13.5		
Welldepth	8 cm		
Welldiameter	4 cm		
Wellvolume	100 ml		