



Hot Bead Sterilizer - 14cm Well / 220 V



This sterilizer provides a safe, effective and convenient method for sterilizing your surgical instruments prior to surgery or tissue culture. Without using potentially dangerous gases, flames or chemicals, you can ensure that your instruments will be free of pathogens and microbial contaminants. The beads take 15-20 minutes to reach 240C to 270C (464F to 518F), and then the sterilizer is ready for use. The temperature is thermostatically maintained, and the unit can be left on all day. Instruments require approximately 20 seconds of contact with the beads for complete sterilization. The sterilizer is safe to use even in laminar flow hoods.

General Info

Product Number	18000-70
Category	Instrument Care > Sterilization
EAN	4251909506148



Find 18000-70 on
finescience.com

Technical Info

Shipping Weight	2450 g	Net Weight	2070 g
Package Size	1 Pc.	Unit of Measure	EA
Minimum Order Quantity	1		
Alloymaterial	Various		
Dimensionstext	20.5 x 12.5 x 13.5		
Voltagepowersource	220 V		
Plugtype	C		
Welldepth	14 cm		
Welldiameter	4 cm		
Wellvolume	175 ml		